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CAUTION!

To work with this microscope, sharp and pointed aids are being used. Please take care that this microscope and its accessories are stored at a place out of reach of children. Let children only work with this microscope under an adult's supervision! Keep packing material (plastic bags etc.) away from children!

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GB Users manual

5 - 8

This manual contains a foldable page at its first and last page. Here are printed all pictures you need for a better understanding of the text.

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GB All parts (fig. 1):

- ① 10x WF Eyepiece
- ③ Barlow lens
- ④ PC-Ocular
- ⑤ Eyepiece holder
- ⑥ Microscope head
- ⑦ Objective-revolver
- ⑧ Microscope stage
- ⑨ Focus wheel
- ⑩ Electric illumination
- ⑪ Electricity supply
- ⑫ Microscope base
- ⓧ Locking screw
- Ⓨ Holding clamps (Slide)

Accessories (fig. 2):

- ① 10x WF Eyepiece
- ② 16x WF Eyepiece
- ③ Barlow lens
- ⑬ Slides, Covering glasses and preparations (5x each) in a plastic box
- ④ PC-Ocular
- ⑬ Ulead Photo Explorer SE Software
- ⑭ AC adapter
- ⑯ Aluminium accessory box
- ⑰ Nonius scale
- ⑱ 5x WF Eyepiece

1. General/Location

Before you set up the microscope, you must choose a suitable location.

First you must make sure that your microscope is on a stable and solid surface.

An electricity supply is required for observation with the electric illuminator.

2. Electric Illumination

It works, by plugging the connection cable into the microscope and the electricity supply. The switch at the rear of the microscope must be switched on.

3. Microscope setup

The microscope's eyepiece (Fig 1, 6) will now be prepared for the first observation.

First, loosen the screw (Fig. 1, X) and rotate the eyepiece into a convenient position.

Begin every observation with the lowest magnification.

Place the microscope's table (Fig. 1, 8) with the focus knob (Fig 1, 9) into the lowest position and rotate the objective revolver (Fig. 1, 7) until it locks on the lowest magnification (4x)

Tip:

First, drive the Microscope's table (Fig. 1, 8) in its lowest position before changing the objective in order to prevent damages.

Important Tip:

According the currently used preparation, high magnifications don't lead in every case to a better picture!

Act very careful here. If you drive the Microscope's table too fast upwards, Objective and slide may touch each other and become damaged!

Insert the 5x eyepiece (18) in the Barlow lens (3).

Take care, that the Barlow lens (2) is inserted completely in the monocular head (3).

4. Observation

After you have set up the microscope with the corresponding illumination, the following principles are important:

Begin each observation with the lowest magnification, so that the centre and position of the object to be viewed is in focus.

The higher the magnification the more light is required for good picture quality.

Begin with a simple observation.

With the mechanical desk (Fig. 3, 17) you can look at your preparation (Fig. 3, 15) in a precise position and to the exact millimeter.

The object is placed between the clamps on the mechanical desk. Move the object, with help of the axis-adjustments, (Fig. 3, X/Y) directly under the objective.

With the built in vernier (Fig. 3, Z) at both axes you can now specifically set and shift the object. It can now viewed with different magnifications.

Look through the eyepiece (1) and turn carefully the focusing wheel (6) until you can see a sharp picture.

Now you can get a higher magnification, while you pull out slowly the Barlow lens (2) of the monocular barrel (3). With nearly entirely pulled out Barlow lens the magnification is raised to 1,6x.

For still higher magnification you can put the 16x eyepiece into the objective revolver and set on higher position (10x / 40x). In this setting the magnification raises also up to 1,6x by pulling out the Barlow lens.

Please notice:

With changing magnification (eyepiece or objective lens changes, pulling out of the Barlow lens) the sharpness of the image must be newly defined by turning the focusing wheel (6).

5. Viewed Object – condition and preparation**Condition**

With an ordinary magnifying glass, we preferably view obscure items like little animals, plants etc. Here light shines on the viewed item, gets reflected and gets through the lens into our eye (Top illumination principle). With a microscope, a so-called Transmission-Type-Microscope, only transparent items can be observed. Light gets through the item, becomes magnified by the objective and the eyepiece and gets in our eye. Many small organisms of the water, plant parts and finest animal components have now from nature these transparent characteristic, other ones must be accordingly prepared. Is it that we make it by means of a pre-treatment or penetration with suitable materials (media) transparent or thus that we cut finest wafers off of them (hand cut, Microtome cut) and these then examine. With these methods will us the following part make familiar.

Manufacture of thin preparation cuts

As already before implemented, as thin cuts as possible are to be manufactured by an object. In order to come to best results, we need something wax or paraffin. If no such material is contained in your microscope set, then you take simply a candle. The wax is given to a pot and warmed up over a flame. The object is dipped now several times into the liquid wax. Let the wax become hard. With a microtome or a knife / scalpel (caution!!!) now finest cuts are cut off by the object coated with wax. These cuts are put on a glass slide and covered with a cover glass.

Manufacture of an own preparation

Put the object which shall be observed on a glass slide and give with a pipette a drop of distilled water on the object (Fig. 7). Set a cover glass (in each well sorted hobby shop available) perpendicularly at the edge of the water drop, so that the water runs along the cover glass edge (Fig. 8). Lower now the cover glass slowly over the water drop.

6. Experiments

If you made yourself familiar with the microscope already, you can accomplish the following experiments and observe the results under your microscope.

6.1 Newspaper print**Objects:**

1. A small piece of paper from a newspaper with parts of a picture and some letters

2. A similar piece of paper from an illustrated magazine
Use your microscope at the lowest magnification and use the preparation of the daily paper. The letters seen are broken out, because the newspaper is printed on raw, inferior paper. Letters of the magazines appear smoother and more complete. The picture of the daily paper consists of many small points, which appear somewhat dirty. The pixels (raster points) of the magazine appear sharply.

6.2 Textile fibers

Items and accessories:

1. Threads of different textiles: Cotton, line, wool, silk, Celanese, nylon etc..
2. Two needles

Each thread is put on a glass slide and frayed with the help of the two needles. The threads are dampened and covered with a cover glass. The microscope is adjusted to a low magnification. Cotton staples are of vegetable origin and look under the microscope like a flat, turned volume. The fibres are thicker and rounder at the edges than in the centre. Cotton staples consist primary of long, collapsed tubes. Linen fibres are also vegetable origin; they are round and run in straight lines direction. The fibres shine like silk and exhibit countless swelling at the fibre pipe. Silk is animal origin and consists of solid fibres of smaller diameter contrary to the hollow vegetable fibres. Each fibre is smooth and even moderate and has the appearance of a small glass rod. Wool fibres are also animal origin; the surface consists of overlapping cases, which appear broken and wavy. If it is possible, compare wool fibres of different weaving mills. Consider thereby the different appearance of the fibres. Experts can determine from it the country of origin of wool. Celanese is like already the name says, artificially manufactured by a long chemical process. All fibres show hard, dark lines on the smooth, shining surface. The fibres ripple themselves/crinkle after drying in the same condition. Observe the thing in common and differences.

6.3 How does bread mushroom develop?

Object: An old piece of bread

The spores of the kind of mushroom, which strike our bread, are to be found everywhere in the atmosphere. Put bread on a slide and squirt carefully some water on it. Moistened bread only, don't wet it. Put the whole into a container with a screw-type cap and to place it into a cabinet, into which only little light breaks in and prevails in a warm temperature. Within a short time the black bread mushroom forms. Regard the bread each day. At the first the mould shows up a white, shining consistence. Take it on a slide to observe it. The material turns out as a complicated thread mass, which forms the mushroom body in its whole. One calls the whole Mycelium. Each thread is a Hyphe. Soon some Rhizoids arise, which embody the mould fungus with bread, in order thereby to water and nutrients for the growth of the Mycelium to receive. In the course of time the Rhizoid colours itself brownish. Vertically over this group Hyphen grow like long slim stacks, which end in a tiny small, white ball. One calls the stack Sporangiohores (carrier of the Sporcap), the ball is a Sporangium or a Sporcapl. Soon these balls accept a black color. Inside sporen present mature. If now the Sporcap breaks open, then it sets the spors free, which step now to air and infect other bread. With the naked eye you can mature the Sporcaps as tiny black marks to recognize. They are scattered on the mould fungus surface and give thus to the kind of mushroom its name. There are however still different kinds of mould fungi. They can be pink, red, blue or green. Manufacture yourselves preparations of all stages of the bread mould.

7. Setup PC-Ocular

Remove the Barlow lens (Fig. 6, 3) with the currently used eyepiece from the tube (Fig. 1, 5) and put the PC-eyepiece (Fig. 5, 4) in place instead (Fig. 5)

8. Installing the software

1. Attach the PC-Ocular to a free USB Port on your computer and turn on your PC.

Please note: It may be possible that the PC-Ocular needs to be attached to the computer AFTER the software installation. In this case boot up your computer without connecting the PC-

Ocular and continue with step 4. To find out which installation method is required by your PC, contact your computer dealer.

2. The Software Serial Number can be found on the CD cover and should be noted as you will need these later during the installation.
3. After boot up, the Hardware Assistant (Fig. A) will pop up on your monitor, indicating that a new device has been attached to your computer and that it will search for a suitable driver. Follow the instructions given by the Hardware Assistant.
4. Insert the software CD into your CD-ROM drive. The software menu (Fig. B) should be loaded up automatically. Click on Install Ulead Photo Explorer 7 SE to commence the installation.

Notice:

Should the CD-Menu not automatically start you should open your Windows-Explorer and select your CD-ROM drive. Open the file <Setup.exe> from the CD and follow the step above.

5. You will now be prompted to select your language of choice. Should your desired language not be available then please select <English> and click on <Next>.
6. The Install Wizard menu appears on your monitor. Click on <Next>.
7. You will now need to accept the <Licence Agreement> by clicking on <Yes> in order to proceed.
8. Please now enter the <User Information>. Enter your name and if applicable your company name and the serial number (found on the CD cover). Now click on <Next> in order to proceed.

Notice:

If the data has not been entered correctly a warning will prompt you to correct the user information before you can continue.

9. The next step will ask you for the destination of where the files should be saved on your harddrive. Click <Next> in order to automatically accept the suggested location.

If you wish to install the software into a different folder you can select it under <Browse>. This folder must be created before the installation to be seen in the menu.

Notice:

The selection of an optional folder can be of significance if your working with a network and would like to install on a drive other than <C:>. In general, we suggest to accept the programs automatic selection for local drives.

10. The installation program will now prompt you under <Create Shortcut Icons>, if various program symbols should be created. Click on all possible options and continue with <Next>.
11. You will now be prompted to create the <Program Folder>. Confirm the name by clicking <Next>.
12. Under <TV System> select <Your country>, and choose the country in which you are living. Continue the installation by pressing <Next>. This step may take several minutes!
13. Under <File Association> select <Associate image file formats with Ulead Instant Viewer> and confirm by pressing <Next>.
14. The software has now been installed. The welcome window of the installation wizard will pop up and the driver installation for the PC-Ocular (ICM532) should now be undertaken by pressing <Next>.

Notice:

It may be possible that complications will occur with older already installed drivers. The computer may show a warning notice informing you to uninstall the old driver before a new one can be installed. You can confirm this step by pressing <OK> and can try working with the old driver. Should the old driver not function correctly then the file <microocular.cat> located on the CD must be installed.

Information on uninstalling and newly installing software can be found in the corresponding section in your Windows-handbook.

15. The window <Select Program Folder> will now appear. Confirm the selection by pressing <Next>.

16. The installation process will be completed.

This will be confirmed with the message <InstallShield Wizard Complete>. Now select <Yes, I want to restart my computer now> and click on <Finish>.

The computer will now be rebooted and the newly installed system data will be updated. This may take up to several minutes.

17. During the reboot process the computer will inform you that the windows driver databank has been updated. Your PC-Ocular is ready for use after the reboot.

Caution:

It may be possible that in some cases the installation has not been properly completed and that the necessary driver for the PC-Ocular can not be found. In this case the driver <microocular.cat> must be manually installed, using the hardware assistant. More detailed information on this topic can be found in your Windows user manual.

9. Working with the PC-Ocular

1. Place a prepared slide onto the microscope stage and properly focus it.

2. Connect the cable of the PC-Ocular to a free USB port of your computer. Remove the eyepiece from the eyepiece holder of your microscope and insert the PC-Ocular instead. Now turn on your computer.

Notice:

The dustcap in front of the PC-Ocular lens must be removed prior to use!

3. After your computer has been properly booted up you can start the ULEAD PhotoExplorer 7 SE software.

4. To view/record streaming video click on the menu option <Video> and then <Capture Video>. You are now in the video mode and can record your observation using the following elements:

- Record
- ▶ Play
- ⏸ Pause
- Stop
- ▶▶ Forward slow / fast
- ◀◀ Rewind slow / fast
- Save as = saving the current recording
- Close = closes the capturing window

Recordings with sound, when selected and the device is attached (PC-Ocular does not support sound recordings).

The recording mode is automatically adjusted to the PC-Ocular and does not require a separate <Setup>.

5. Still pictures or picture sequences can be recorded by entering the menu <Video> and clicking on <Capture still image>.

- Live View = shows the current observation
- Save as = save current recording
- Capture Now = creates a single capture, if  is selected
creates a sequence of images (every 5 seconds),
if  is selected
- Close = closes the capturing window

The recording mode is automatically adjusted to the PC-Ocular and does not require a separate <Setup>.

A more detailed instruction manual of the Ulead PhotoExplorer 7 SE in various different languages is located on the software CD as a <pdf> document. Should your language not be available on the CD contact Ulead directly to see if your language is available.

10. Care and maintenance

The microscope is a high quality optical appliance. Therefore you should make sure that dust or moisture do not come into contact with your microscope.

Avoid putting fingerprints on any optical surfaces. If dust or dirt should nevertheless have gotten onto your microscope, or the accessories, you should first remove this with a soft brush. Then clean the soiled area with a soft, lint-free cloth. To remove fingerprints from the optical surfaces it is best to use a lint-free, soft cloth, to which you have applied some alcohol.

After usage, you should stow the microscope and the accessories in their correct containers again.

Remember:

A well maintained microscope will keep its optical quality for years and thus maintain its value.

11. Troubleshooting

Error	Solution
No picture visible	switch light on / adjust focus
Picture flickers while viewing with PC eyepiece	if necessary, adjust resolution of the video graphics board
Software installation reports "not XP approved"	confirm by clicking "OK"

12. Specifications

System requirements for PC eyepiece
 Intel Pentium with 166 MHz or equiv. / higher
 Operating System Windows 98 SE / ME / 2000 / XP
 Min. 64 MB RAM
 Min. 100 MB free hard disk space

Magnification table

<u>Eyepiece</u>	<u>Objective</u>	<u>Magnification</u>	<u>(with Barlow lens)</u>
10x	4x	40x	64x
10x	10x	100x	160x
10x	40x	400x	640x
16x	4x	64x	102x
16x	10x	160x	256x
16x	40x	640x	1024x

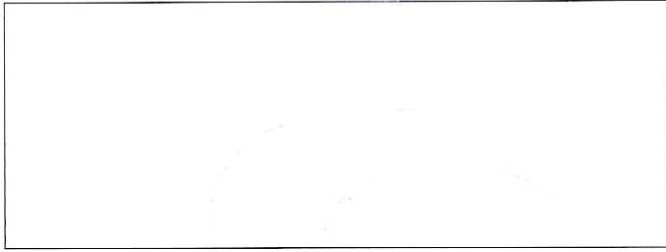
13. Warranty

The period of warranty is two years, beginning on the day of purchase. Please keep the cash receipt as evidence of purchase. Microscopes which become defective during the warranty period can be returned to the dealer where the microscope was bought. The repaired microscopes or a new pair will then be returned to you, free of charge. In the case of defects which occur after the end of the warranty period, the microscopes can also be returned. However, repairs which become necessary after the end of the warranty period will be subject to a service fee.

Important:

Make sure to return the microscopes carefully packed in the original packaging in order to prevent transport damage. Please also enclose the cash receipt (or a copy). This guarantee does not imply any restriction of your statutory rights.

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